

ECT Step by Step

Mycotoxin

8/08/08

Mycotoxin Check Samples

For New Test
Select FGIS Mycotoxin from
Dropdown next to Create New Test

To finish an Open test use Search
or enter the Test Reference No.

You are here: ECT Menu 1.0.7

Checktest Home

Create New Checktest

[Search For Existing Checktest](#)

Search By Test Reference No.

[Equipment](#)

[Reports](#)

[E-mail Distribution Lists](#)

[Administration](#)

- FGIS-904 Laboratory Scale Test
- FGIS-923 Moisture Meter Test
- FGIS-924 Barley Pearler
- FGIS-924 Sieve Test
- FGIS-925 Rice Checktest Form
- FGIS-927 Test Weight Check Test
- FGIS-928 Dockage Check Test
- FGIS-936 Sampler Condition Report
- FGIS-965 Scale Test Report (Hopper)
- FGIS-965-1 Scale Test Report (Railroad Track)
- FGIS-965-2 Scale Test Report (Vehicle)
- FGIS-Falling Number
- FGIS-Mycotoxin**
- FGIS-NIRT Daily Barley
- FGIS-NIRT Daily Corn
- FGIS-NIRT Daily Soybean
- FGIS-NIRT Daily Wheat

Blank General Page

Steps to Complete:

1. Enter Service Point or use Lookup
2. Click on Save
3. Click on Next

You are here: ECT Menu > FGIS-Mycotoxin General Information 1.0.7

**FGIS-Mycotoxin
General Information**

General Standard Test Summary

Test Reference Number:

Service Point Lookup Service Point

FD OA LOCATION

Save Cancel Void Next

Completed General Page

General page must be saved to move to Next page.

You are here: ECT Menu > FGIS-Mycotoxin General Information 1.0.7

**FGIS-Mycotoxin
General Information**

General | Standard | Test | Summary

Test Reference Number: 14244

Data has been saved successfully

Service Point Lookup Service Point

FO	OA	LOCATION
FGIS - Wichita Field Office		KANSAS CITY, MO

Save | Cancel | Void | Next

Last change made on 7/26/2008 10:17:05 PM by user Mary Vick.

Blank Standard Page

1. Select Test Type
2. Select Quarter and Year
3. Enter Sample ID values and Targets.
4. Enter Results By and Date
5. Enter Remarks as needed
6. Click on Save
7. Review
8. Click on Submit or Submit/Create Batch

**FGIS-Mycotoxin
Standard Information**

Test Reference Number: 14244

Test Type
For Qtr Year

Column	Sample ID	Target
1	<input type="text"/>	<input type="text"/>
2	<input type="text"/>	<input type="text"/>
3	<input type="text"/>	<input type="text"/>
4	<input type="text"/>	<input type="text"/>

Results By Date

Remarks

Completed Standard Page

You are here: ECT Menu > FGIS-Mycotoxin Standard Information

1.0.7

FGIS-Mycotoxin Standard Information

General Standard Test Summary

Test Reference Number: 14244

Data has been successfully saved

Test Type Don

For Qtr 2nd Year 2008

Column	Sample ID	Target
1	100	0.0
2	200	25.0
3	300	25.0
4	400	0.0

Results By 94110 Date 7/25/2008

Remarks

Previous

Save

Cancel

Void

Submit

Submit/Create Batch

Next

Batch page

Enter Service Point number and select number of samples for that location.

Submit to complete batch process.

You are here: ECT Menu > Create Checktest Batch 1.0.7

Create Checktest Batch

Test Reference Number: 14244

Service Point		Number of Samples
Service Point <input type="text" value="244970"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		
Service Point <input type="text" value="244300"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		
Service Point <input type="text" value="263620"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		
Service Point <input type="text" value="262830"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		
Service Point <input type="text" value="155400"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		
Service Point <input type="text" value="121240"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		
Service Point <input type="text" value="128530"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		
Service Point <input type="text" value="247070"/> <input type="button" value="Lookup Service Point"/>		<input type="text" value="1"/> ▼
FO OA LOCATION		

Blank Test Page

1. Select Test Kit.
2. Enter Actual Test values.
3. Kit Lot # required or None if unknown.
4. Enter Results By and Date
5. Click on Save
6. Review
7. Click on Submit

You are here: ECT Menu > FGIS-Mycotoxin Test Information

1.0.7

FGIS-Mycoing Number Test Information

General Standard Test Summary

Test Reference Number: 14244

Equipment Type Don

Quarter & Calendar Year 2nd Quarter 2008

Test Kit	Sample ID				Kit Lot #
	100	200	300	400	
<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>
<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>
<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>
<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>
<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>	<input type="text"/>

Results By Date

Remarks

Previous

Save

Cancel

Void

Submit

Next

Completed Test Page

On submit screen moves automatically to Summary Page

FGIS-Mycoing Number Test Information

General Standard Test Summary

Test Reference Number: 14244

Data has been successfully saved

Equipment Type Don

Quarter & Calendar Year 2nd Quarter 2008

Test Kit	Sample ID				Kit Lot #
	100	200	300	400	
Don Fluoroquant	0	17	20	0	ab39438023
Don Ridascreen Fast	0	10	30	1	z3bd54234

Results By 94110 Date 7/28/08

Remarks

Previous Save Cancel Void Submit Next

Summary Screen

Results of test kits and Target values shown

Click on Cancel to go back to Home page or use breadcrumb menu at the top

You are here: ECT Menu > FGIS-Mycotoxin Summary Information

1.0.7

FGIS-Mycotoxin Summary Information

General Standard Test Summary

Test Reference Number: 14244

Test data submitted successfully. Checktest is now closed.

Equipment Type Don

Quarter & Calendar Year 2nd Quarter 2008

Test Kit	Sample ID				Kit Lot #
	100	200	300	400	
Don Fluoroquant	0.0	17.0	20.0	0.0	ab39438023
Don Ridascreen Fast	0.0	10.0	30.0	1.0	z3bd54234
Target	0.0	25.0	25.0	0.0	

Date Tested 7/28/2008

Remarks

Previous Save Cancel Void Submit


Remarks

Enter any additional information in remarks and hit Save.

All remarks are saved the same way. It shows who and when each remarks entry is made.

Remarks never locked.

Remarks



7/9/2008 1:11:44 PM : Mary Vick : Kettle tested with apparatus FM#1

Previous Save Cancel Void

Notes:

A large, empty rounded rectangular box with a dark blue border, intended for taking notes. The box is centered on the slide and occupies most of the lower two-thirds of the page.